




Cheminert® Flangeless Fittings with collapsible ferrule

Cheminert flangeless fittings now include our new patented* collapsible ferrule design. This innovative design utilizes a one-piece ferrule engineered to collapse as it is tightened. The collapse takes place in two very narrow areas, and results in a very effective seal with virtually no distortion of the tubing ID and no dead volume.

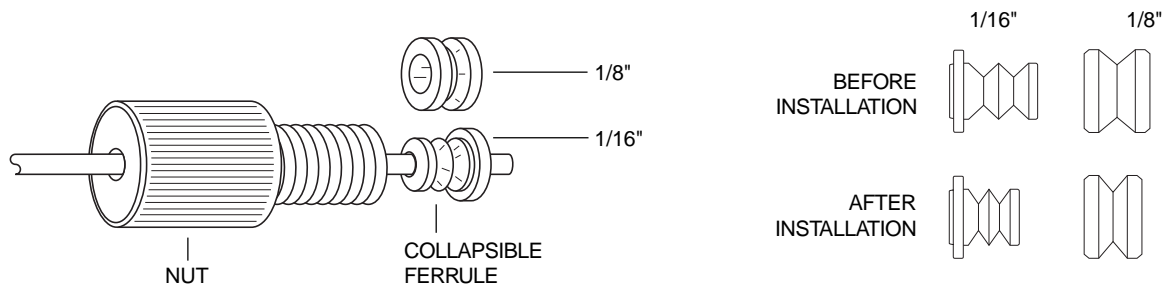

Instructions

1. Use a sharp razor blade (or polymer tubing cutter) to cut the tubing, making sure that the cut is square and that the tubing is not crushed.
2. Slide the nut onto the tubing, followed by the ferrule. With 1/16" ferrules, the small end goes toward the nut. For 1/8" ferrules, the orientation doesn't matter.
3. Screw the assembly part way into a 1/4-28 female fitting (or a Cheminert setting tool, p/n CST).
4. Make sure the tubing is pushed all the way into the female fitting, then screw in the nut until you feel the increase in resistance which signifies that the nut has contacted the ferrule. Finger-tighten it 3/4 turn past that point.

NOTE: These instructions are valid for the standard CTFE collapsible ferrules (p/n CFL-CB1KF or CFL-CB2KF) as well as the optional PEEK ferrules (p/n CFL-CB1PK or CFL-CB2PK), but the PEEK ferrules require a little more force to apply the 3/4 turn. Either ferrule used with PEEK tubing may require one full turn, but *tightening in excess of one turn is not advised*. We strongly recommend the use of PEEK ferrules with PEEK tubing.

Caution:

Extreme overtightening can result in loss of sealing capacity.

Retightening a previously made up fitting

Screw the nut into the fitting detail approximately 1/4 turn after the nut begins to compress the ferrule.

*Patent No. 6,575,501

North America, South America, and Australia/Oceania contact:

VICI® Valco Instruments Co. Inc.
 P.O. Box 55603
 Houston, TX 77255
 Sales: (800) 367-8424
 Tech: (713) 688-9345
 Fax: (713) 688-8106 valco@vici.com

Europe, Asia, and Africa contact:

VICI® VICI AG International
 Parkstrasse 2
 CH-6214 Schenkon
 Switzerland
 Phone: +41 41 925 6200
 Fax: +41 41 925 6201 info@vici.ch

Cheminert® and VICI® are registered trademarks of Valco Instruments Co. Inc. and VICI AG

TN-505a Rev 10/03