




## Universal Actuator Tutorial


### Part 1: Downloading the USB Drivers

Many users may be proficient at downloading and installing drivers in a Windows environment. If that is your situation you may wish to skip parts 1 and 2 of the tutorial and go directly to part 3, "Verifying Actuator Communication."

1. Open your browser and navigate to [www.vici.com](http://www.vici.com).
2. Move your mouse over **PRODUCTS** in the top navigation bar to open the product menu.

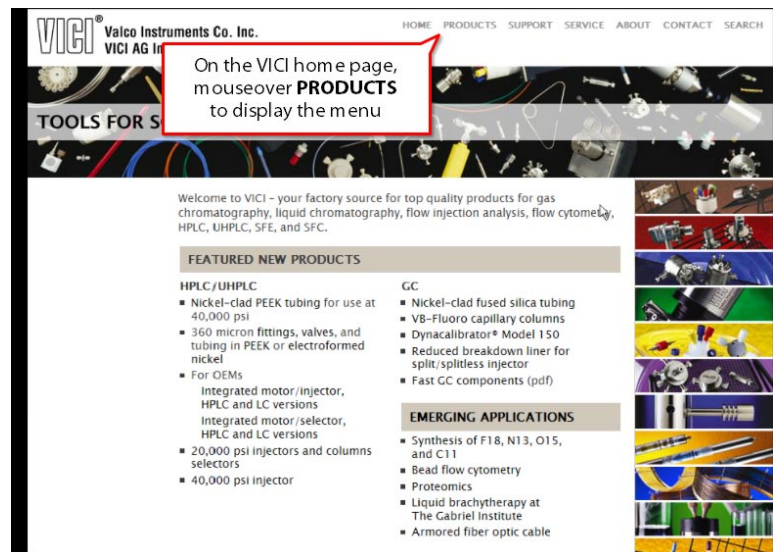

3. Under **ACTUATORS**, select **Universal**.

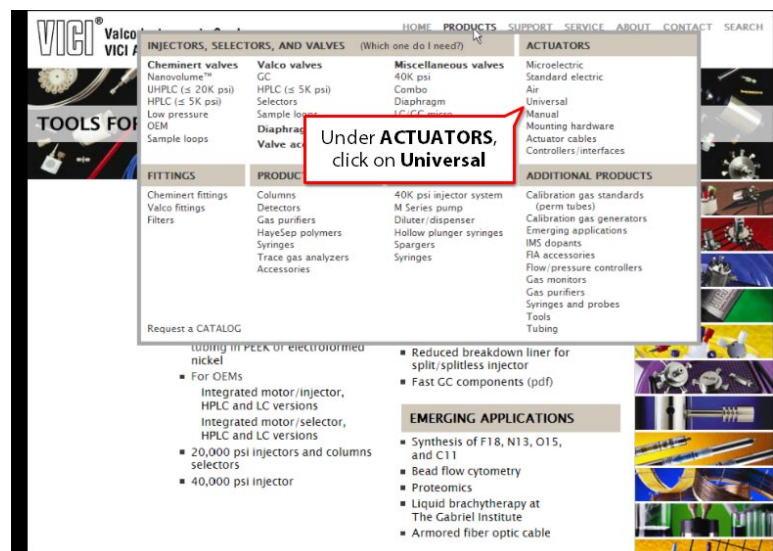

- On the universal actuator page, click on the **USB drivers** link under **MORE INFORMATION**.

**Under MORE INFORMATION, click on USB drivers**

**MORE INFORMATION**

- Which actuator model do I need for the valve I have?
- User manual
- USB drivers (Microsoft WHQL certified)
- Product brochure
- Declaration of CE compliance

**MOUNTING HARDWARE**

Mounting hardware must be ordered separately. Please contact our sales or technical staff to determine the correct hardware needed.

Model	Interface	Product No.	Price
EUH High speed	none	EUH	US customers <a href="#">log in</a> for prices.
	RS-232	EUHA	All others <a href="#">click here</a> for purchasing information.
	USB	EUHB	
	BCD	EUHC	

- When prompted, choose to save the file, rather than opening it.
- Then click **OK**.

**Make sure the Save File radio button is selected**

**Then click OK**

**MORE INFORMATION**

- Which actuator model do I need for the valve I have?
- User manual
- USB drivers (Microsoft WHQL certified)
- Product brochure
- Declaration of CE compliance

**MOUNTING HARDWARE**

Mounting hardware must be ordered separately. Please contact our sales or technical staff to determine the correct hardware needed.

Model	Interface	Product No.	Price
EUH High speed	none	EUH	US customers <a href="#">log in</a> for prices.
	RS-232	EUHA	All others <a href="#">click here</a> for purchasing information.
	USB	EUHB	
	BCD	EUHC	

- Select an appropriate download location. Often that will be Downloads or My Downloads – just make sure to remember where it's getting saved.
- Click **Save**.

**Check to see that the location is appropriate**

**Then click SAVE**

**MORE INFORMATION**

- Which actuator model do I need for the valve I have?
- User manual
- USB drivers (Microsoft WHQL certified)
- Product brochure
- Declaration of CE compliance

**MOUNTING HARDWARE**

Mounting hardware must be ordered separately. Please contact our sales or technical staff to determine the correct hardware needed.

Model	Interface	Product No.	Price
EUH High speed	none	EUH	US customers <a href="#">log in</a> for prices.
	RS-232	EUHA	All others <a href="#">click here</a> for purchasing information.
	USB	EUHB	
	BCD	EUHC	