




## Universal Actuator Tutorial

### Part 2: Installing the USB Drivers


Many users may be proficient at downloading and installing drivers in a Windows environment. If that is your situation you may wish to skip parts 1 and 2 of the tutorial and go directly to part 3, "Verifying Actuator Communication."




1. Click **start** to begin. When the options expand, click on **My Computer**.

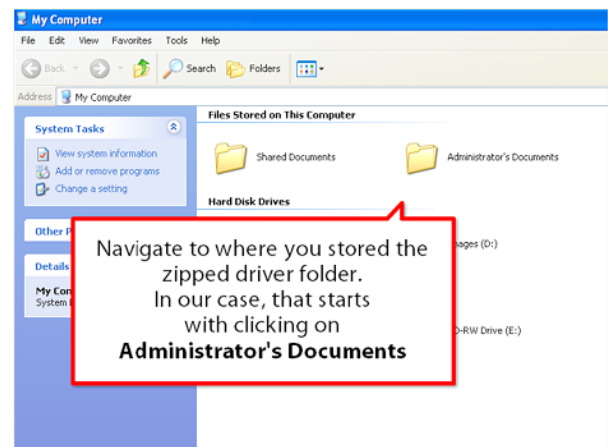


2. Navigate to the location you choose to save the compressed driver folder in part 1 of the tutorial.

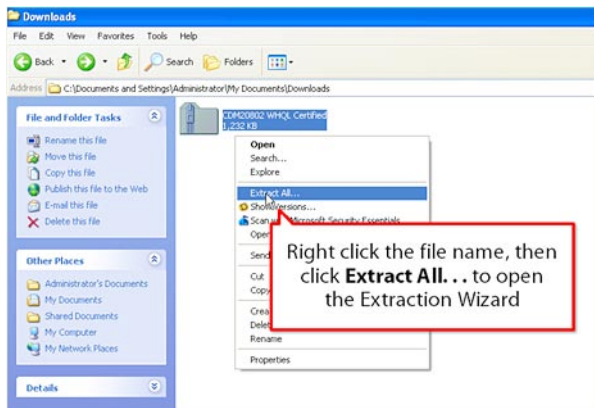


3. Right click on the compressed folder, then select **Extract All** to open the Extraction Wizard. When the wizard opens, click **Next**.

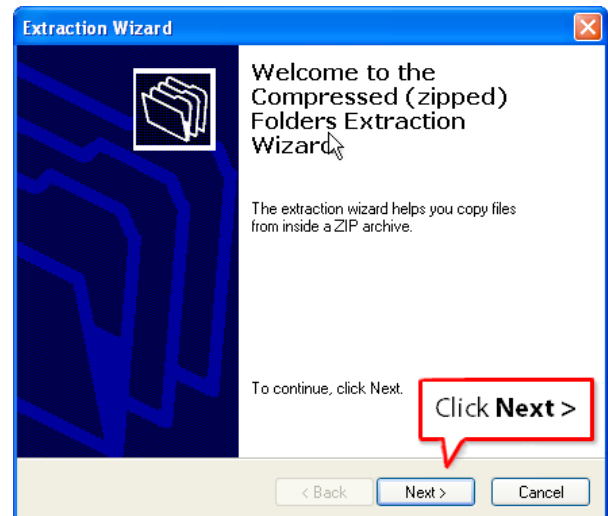


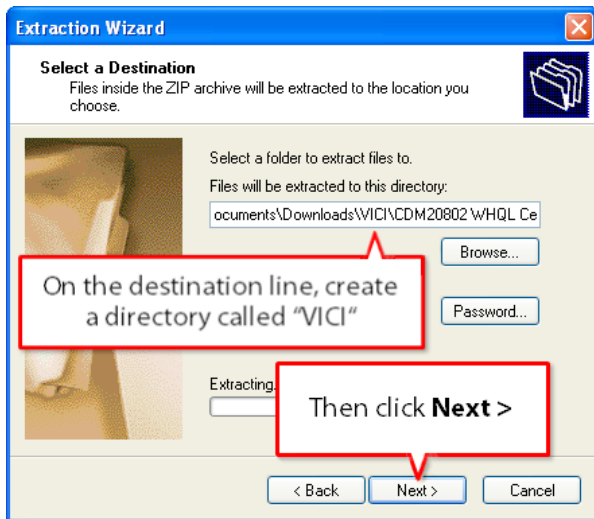


4. Type "VICI" in front of the name of the compressed folder on the destination line. When you click **Next**, the files will be unzipped to the newly-created VICI directory.

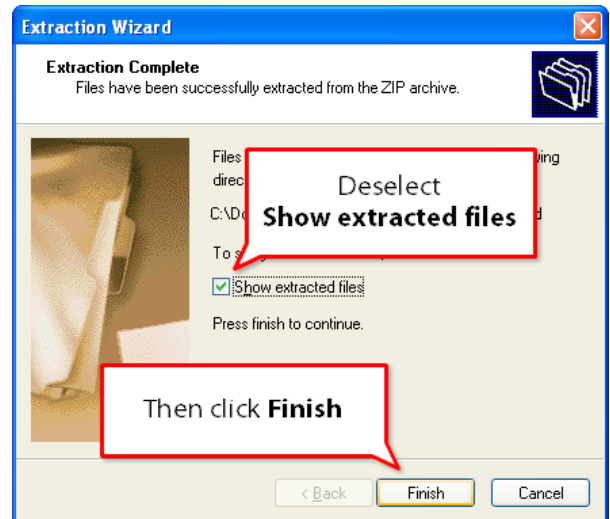


5. Since there is no need to view the extracted files, deselect **Show extracted files** and click **Finish**.

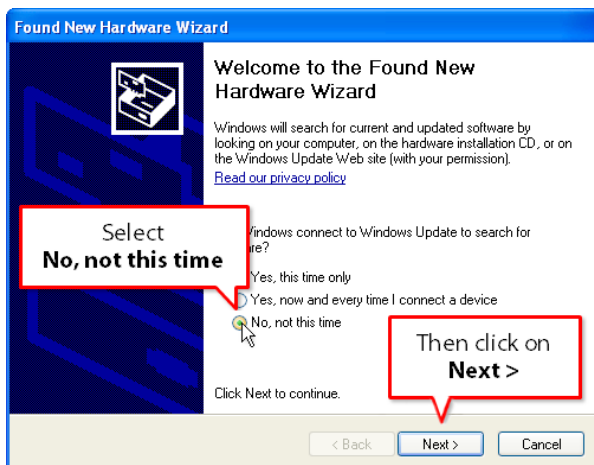


6. Connect the power cable to the universal actuator and plug it to an appropriate electrical outlet.  
7. Plug the USB cable into the actuator and into your computer. After a few moments, the screen above will appear. Select **No, not this time**, then click **Next**.

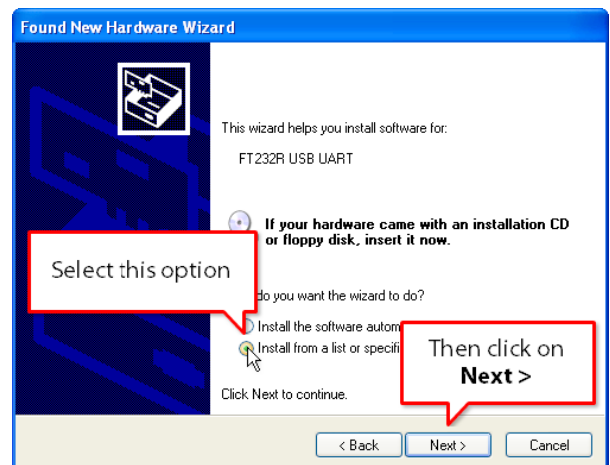


8. Select **Install from a list or specific location**, then click **Next**.

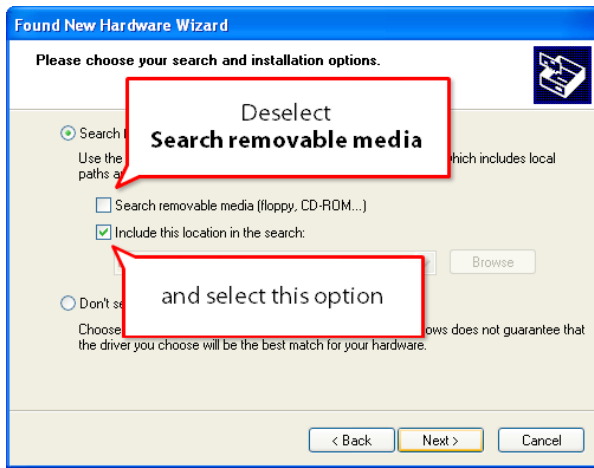


9. Deselect **Search removable media** and select **Include this location in the search.**

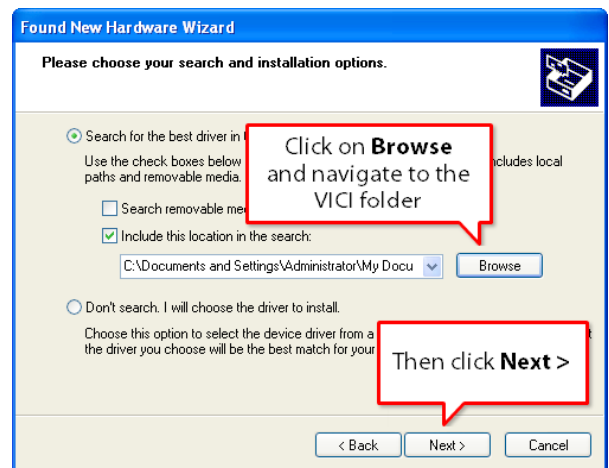

10. Click on **Browse** to navigate to the VICI folder containing the uncompressed driver files. Then click **Next**.

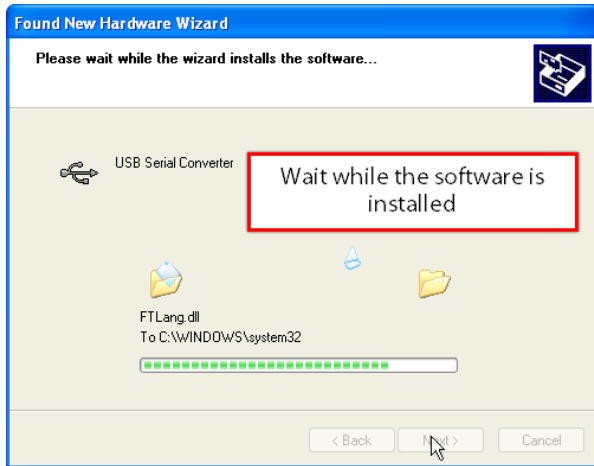

11. A series of progress screens indicate that the installation is underway.

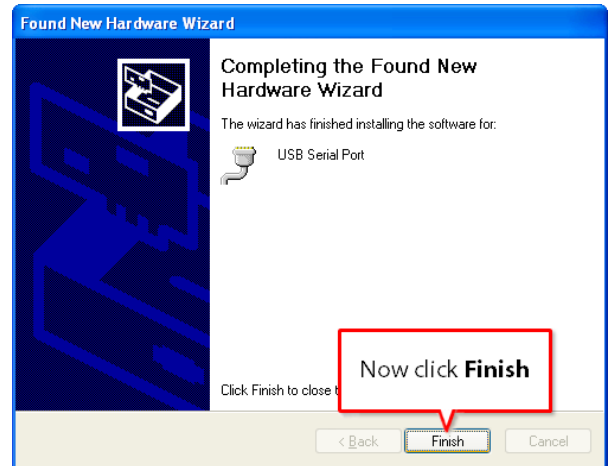

12. When the installation is complete, click **Finish**, and you're done!